

Shut Down and clean up procedures

When you are done imaging, remove your sample from the microscope and place it on a kimwipe or other paper towel if it has oil or water.

If you used an oil objective, please refer to the [instructions](#) of Dr. Dragavon to clean it.

Shut down the system according to the shut-down procedure taped on the wall:

- A. Save your beautiful images to your shared drive.
- B. Exit Nikon Elements and disconnect from your shared drive.
- C. Remove your sample from the stage and clean up the microscope (remove oil from the objective, shutdown the environmental chamber...). Manually lower the objective to a position below 1000 on the LCD screen.
- D. Open the reservation calendar. If there is a user within 2 hours of your session, proceed to F. If the user is more than 2 hours after your session, proceed to Step E.
- E. Turn the key on the laser box to the LEFT (3).
- F. Turn off the two power strips (1 and 2).
- G. Publish an awesome article.

Revision #1

Created 9 October 2023 17:51:11 by Evolene Premillieu

Updated 9 October 2023 17:53:56 by Evolene Premillieu